Faustovirus Capping Enzyme (S17)

Code No. 2480A **Size:** **500 U**
Conc.: **25 U/ μ l**

Supplied Reagents:

10X Capping Buffer 2 **100 μ l**
S-adenosylmethionine (SAM) (32 mM) **100 μ l**
GTP (10 mM) **50 μ l**

Description:

Faustovirus Capping Enzyme (S17) is a single-subunit enzyme derived from the Faustovirus S17 strain, which adds a 7-methylguanylate structure (Cap 0) to the 5' end of RNA (5'-triphosphate RNA) in a similar manner to the action of Vaccinia virus capping enzyme. In eukaryotes, cap structures play crucial roles in stabilizing, nuclear exporting, and translating mRNAs. By using this enzyme in conjunction with the provided reagents (10X Capping Buffer 2, S-adenosylmethionine (SAM), and GTP), the efficient production of intact Cap-0 RNA can be achieved *in vitro*. Furthermore, when combined with mRNA Cap 2'-O-Methyltransferase (Cat. #2470A/B), this enzyme facilitates the production of intact Cap-1 RNA, which features a methyl group at the 2'-O position of the first nucleotide of Cap-0 RNA, in a single-step reaction. Moreover, compared to a Vaccinia virus capping enzyme, this enzyme exhibits a wider temperature range for its enzymatic activity (30 - 50°C) and is more versatile in working with various targets. These characteristics enable the robust preparation of mRNA for use in research and development in the field of RNA therapeutics, such as vaccines.

Storage: -20°C

Source:

Escherichia coli carrying a plasmid containing the gene for Faustovirus capping enzyme (S17)

Property:

Molecular mass: approx. 100 kDa

Unit Definition:

One unit is defined as the amount of enzyme that converts 5 pmol of a 100 nt transcript to Cap-0 RNA in 30 minutes at 37°C.

Reaction Mixture for Unit Definition:

1X	Capping Buffer 2
25 μ M	GTP
20 μ Ci/ml	[3H] SAM
2.5 μ g/50 μ l	100 nt RNA

Quality Control Data:

Please see the Certificate of Analysis (CoA) for each lot. You can download the CoA on Takara Bio website.

Applications:

1. Preparation of Cap-0 mRNA
2. Preparation of Cap-1 mRNA in conjunction with mRNA Cap 2'-O-Methyltransferase

Precautions for Use:

1. Do not mix the enzyme vigorously.
2. RNA is less produced or fragmented when reagents, tubes, or pipette tips are contaminated with RNase. Precautions should be taken during experiments to avoid RNase contamination, such as wearing disposable gloves and using tubes and tips exclusively for RNA experiments.

Protocols:

1. Preparation of Cap-0 RNA from uncapped RNA (5'-triphosphate RNA)

RNA transcript (50 μ g)*1	38 μ l
10X Capping Buffer 2	5 μ l
GTP (10 mM)	2.5 μ l
SAM (2 mM, dilute 32 mM stock to 2 mM)*2	2.5 μ l
Faustovirus Capping Enzyme (S17) (25 U/ μ l)*3	2 μ l
Total	50 μl*4

Incubate at 37°C for 30 minutes.*3

2. Preparation of Cap-1 RNA from uncapped RNA (5'-triphosphate RNA)

RNA transcript (50 μ g)*1	34 μ l
10X Capping Buffer 2*5	5 μ l
GTP (10 mM)	2.5 μ l
SAM (4 mM, dilute 32 mM stock to 4 mM)*2	2.5 μ l
Faustovirus Capping Enzyme (S17) (25 U/ μ l)*3	2 μ l
mRNA Cap 2'-O-Methyltransferase (50 U/ μ l)*3	4 μ l
Total	50 μl*4

Incubate at 37°C for 1 hour.*3

- *1 The heat denaturation of the RNA (5'-triphosphate RNA) prior to the reaction may increase the capping efficiency by reducing the secondary structure at the 5' end of the transcript.
 - (1) Take 50 μ g of the RNA and adjust the volume to 38 μ l (Cap-0 RNA prep) or 34 μ l (Cap-1 RNA prep) with RNase-free Water.
 - (2) Heat-denature the RNA at 65°C for 5 - 10 minutes, then immediately place it on ice for 5 minutes.
- *2 SAM is unstable. Dilute the necessary amount of the 32 mM stock solution with RNase-free Water prior to the reaction, and keep it on ice until use.
- *3 If the capping efficiency of RNA is low, increase the amount of enzymes used and/or extend the incubation time.
- *4 The reaction can be scaled up as needed.
- *5 Use this buffer provided for Faustovirus Capping Enzyme (S17) instead of the buffer provided with mRNA Cap 2'-O-Methyltransferase (Cat. #2470A/B).

Related Products:

mRNA Cap 2'-O-Methyltransferase (Cat. #2470A/B)

RNase-free Water (Cat. #9012)

Takara IVTpro™ mRNA Synthesis System (Cat. #6141)

Cloning Kit for mRNA Template (Cat. #6143)

Takara IVTpro™ T7 mRNA Synthesis Kit (Cat. #6144)

Template Vector (*BspQ I*) for T7 mRNA Synthesis (Cat. #6146)

BspQ I (Cat. #1227A/B)

T7 RNA Polymerase ver.2.0 (Cat. #2541A)

PrimeCap™ T7 RNA Polymerase (low dsRNA) (Cat. #2560A)

Recombinant RNase Inhibitor ver.2.0 (Cat. #2315A/B)

Pyrophosphatase (inorganic) (Cat. #2450A/B)

IVTpro and PrimeCap are trademarks of Takara Bio Inc.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us by our website at www.takarabio.com.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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2. Uncapped RNA (5'-triphosphate RNA) から Cap-1 RNA の調製

RNA transcript (50 μ g) *1	34 μ l
10×Capping Buffer 2 *5	5 μ l
GTP (10 mM)	2.5 μ l
SAM (4 mM, dilute 32 mM stock to 4 mM) *2	2.5 μ l
Faustovirus Capping Enzyme (S17) (25 U/ μ l) *3	2 μ l
mRNA Cap 2'-O-Methyltransferase (50 U/ μ l) *3	4 μ l
Total	50 μ l *4

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