α -1,3/4-L-Fucosidase

Code No. 4453

Size: 100 μ U/100 μ l

Systematic name: α -L-Fucoside fucohydrolase

Enzyme code: 3.2.1.51

Source: *Streptomyces* sp. 142

Form:
Solution in 50 mM potassium phosphate buffer, pH 6.0, containing 0.1 M NaCl, 0.02% NaN₃ and 0.1% Brij-58.

Storage: 4°C

Reaction:
This enzyme catalyses the hydrolytic cleavage of α -fucoside bond and the L-fucose is released.

Substrate specificity:

- This enzyme specifically cleaves α 1, 3- and α 1, 4-fucoside bond.
- This enzyme also releases L-fucose from α 1-acid glycoprotein.

Properties:

Molecular weight: 40,000 (Gel filtration method)

Michaelis constant:

- K_m=10.1 μ M (PA-Lacto-*N*-fucopentaose II: Table 1, Substrate 1)
- K_m=1.08 μ M (PA-Lacto-*N*-fucopentaose III: Table 1, Substrate 2)

Optimum pH: pH 6.0 (Table 1, Substrate 2)

Range of pH stability: pH 4.5 - 7.0 (37°C, 60 minutes)

Stabilizer: 0.1% NP-40, 0.1% Brij-58, and 0.02% BSA

Definition of activity:
One unit is the amount of enzyme required to hydrolyze 1 μ mol of pyridylamino lacto-*N*-fucopentaose III within 1 minute at 37°C at pH 6.0.

Purity:

<Protease Contamination>

No protease activity is detected after incubation of 10 μ U α -1,3/4-L-Fucosidase with 4 nmol oxidized insulin B chain for 16 hours at 37°C, in 20 μ l of 150 mM phosphate buffer, pH 7.5.

<Exoglucosidase Contamination>

No detectable α -galactosidase, β -galactosidase, β -*N*-acetylglucosaminidase, α -mannosidase, and β -xylosidase activities are detected when 20 μ U α -1,3/4-L-Fucosidase is incubated with 100 nmol of a *p*-nitrophenyl glycoside for 16 hours at 37°C, in 0.1 ml of 300 mM potassium phosphate buffer, pH 6.0.

References:

- 1) Sano M, Hayakawa K, and Kato I. *J Biol Chem.* (1992) **267**: 1522-1527.
- 2) Kondo A, Suzuki J, Kuraya N, Hase S, Kato I, and Ikenaka T. *Agric Biol Chem.* (1990) **54**: 2169-2170.
- 3) Suzuki J, Kondo A, Kato I, Hase S, and Ikenaka T. *Agric Biol Chem.* (1991) **55**: 283-284.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com. Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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PA: Pyridylamino (group) PA-Sugar chain was derivatized according to the method of Kondo *et al.*

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* : 4 M trifluoroacetic acid 100°C、3 hr

α -1,3/4-L-Fucosidase

Application 1 : Enzymatic cleavage of PA-Sugar chain with α -1,3/4-L-Fucosidase

The reactions were performed using PA-sugar chains in Table 1 as substrates.

The reaction mixture was as follows :

10 μ M PA-Sugar chain	2 μ l
500 mM Potassium phosphate buffer, pH 6.0	2 μ l
Enzyme solution	X μ l
Sterile purified water	6 - X μ l
Total	10 μ l

The reaction mixture was incubated at 37°C for 30 min or 60 min, then 40 μ l of 1% trifluoroacetic acid was added to reaction mixture to stop the reaction. The reaction mixture was analyzed by HPLC.

Result :

Substrate (Number of Table 1)	Amount of enzyme (μ U)	Cleaved product	Rate of hydrolysis (%)	
			30 min	60 min
1	1	Gal β 1-3GlcNAc β 1-3Gal β 1-4Glc-PA	100	100
2	1	Gal β 1-4GlcNAc β 1-3Gal β 1-4Glc-PA	95	100
3	5	Fuc α 1-2Gal β 1-3GlcNAc β 1-3Gal β 1-4Glc-PA	77	100
6	2	 Gal β 1-4GlcNAc β 1-2Man α 1 Gal β 1-4GlcNAc β 1-4Man α 1 Gal β 1-4GlcNAc β 1-2Man α 1 Man β 1-4GlcNAc β 1-4GlcNAc-PA	91	100
8	5	Fuc α 1-2Gal β 1-3GlcNAc β 1-3Gal β 1-4Glc-PA	0	0

Application 2 : Enzymatic cleavage of glycoprotein with α -1,3/4-L-Fucosidase

α 1-acid glycoprotein (α 1-AGP) is the serum glycoprotein, M.W. 44,000 and it contains approx. 45% of sugar in the molecule. Its sugar chains are linked to five intramolecular Asn residues by *N*-glycoside bond and its structures are *N*-Acetylglucosamin type biantennary, triantennary, and tetraantennary sugar chains.

Since a fucose is linked to the nonreducing terminal GlcNAc of triantennary or tetraantennary unit by α -1,3 bond, α -1,3/4-L-Fucosidase was used to release the fucose.

50 pmol of α 1-AGF and 25 μ U of α -1,3/4-L-Fucosidase was incubated at 37°C for 20 hours in 10 μ l of 100 mM potassium phosphate buffer, pH 6.0, containing 2 M ammonium sulfate, then 1% trifluoroacetic acid was added to reaction mixture to stop the reaction.

20 μ l of reaction solution was dried, then the released fucose was pyridylaminated by using PALSTATION. PA-fucose was determined by HPLC using the anion exchange column (PALPAK Type A).³⁾

Result :

The activity of this enzyme for glycoprotein was increased by adding ammonium sulfate remarkably.

Cleavage Condition		Amount of released L-Fucose (%)
Acid Hydrolysis *		100
Enzymatic Cleavage	(NH ₄) ₂ SO ₄ , (-)	11
	(NH ₄) ₂ SO ₄ , (+)	101

* 4 M trifluoroacetic acid 100°C, 3hr