LA PCR™ Genome DNA Set

Code No. 9060

Size: ea. 20 reactions

Description:

This set includes highly purified, high-molecular DNA originated from human genome and *E. coli* genome to be used as the template for LA PCR. Also included are control primers to amplify 17.5 kb fragment using human genome DNA as the template, and 20 kb using *E. coli* genome DNA.

Components:

human genome DNA	10 μ g (500 ng/ μ l)
<i>E. coli</i> genome DNA	2 μ g (100 ng/ μ l)
human-1 primer	200 pmol (10 pmol/ μ l)
human-2 primer	200 pmol (10 pmol/ μ l)
<i>E. coli</i> -1 primer	200 pmol (10 pmol/ μ l)
<i>E. coli</i> -2 primer	200 pmol (10 pmol/ μ l)

Storage:

DNA : 4°C
(It can be stored at -20°C . Avoid freeze-thaw cycle.)

Primer : -20°C

Form:

DNA :	human :	10 mM Tris-HCl, pH 8.0
		0.1 mM EDTA
	<i>E. coli</i> :	10 mM Tris-HCl, pH 8.0
		1 mM EDTA
Primer:	ea.	10 mM Tris-HCl, pH 8.0
		0.1 mM EDTA

Origin:

human genome DNA : HL60
(Human promyelocytic leukemia cell)

E. coli genome DNA : JM109

Purity:

human genome DNA : $A_{260}/A_{280} > 1.8$

E. coli genome DNA : $A_{260}/A_{280} > 1.8$

Primer :

human-1 :	5'-ACATGATTAGCAAAGGGCCTAGCTTGGACTCAGA-3'
human-2 :	5'-TGCACCTGCTCTGTGATTATGACTATCCCACAGTC-3'
<i>E. coli</i> -1 :	5'-GGTGGCGATGCAAATGCAATCTTCGTTGCCCAAC-3'
<i>E. coli</i> -2 :	5'-TGCAAATACTTCTGCGCAATGCGGCATTGAAGT-3'

General reaction mixture for PCR :

Takara LA Taq® (5 U/ μ l)	0.5 μ l
10X LA PCR Buffer II (Mg ²⁺ free)	5 μ l
25 mM MgCl ₂	5 μ l
dNTP Mixture (2.5 mM each)	8 μ l
human genome DNA (500 ng/ μ l) or <i>E. coli</i> genome DNA (100 ng/ μ l)	1 μ l
Primer 1 *	1 μ l or 0.2 μ M
Primer 2 *	1 μ l or 0.2 μ M
Sterile Purified water	up to 50 μ l

* In case using control primer set, use 1 μ l each.

Conditions of thermal cycling (In case of control primer set) :

94°C	1 min	} 30 cycles
98°C	10 sec	
68°C	15 min	
72°C	10 min	

Reference :

Silhavy T J, *et al.* Experiments with Gene Fusions.
Cold Spring Harbor Lab. Cold spring Harbor N.Y. (1984) 137-139.

Takara LA Taq is a registered trademark of Takara Bio Inc.
LA PCR is a trademark of Takara Bio Inc.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com. Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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